

! micrux  
TECHNOLOGIES

---




Thick-film Carbon  
Single-Electrodes



**Thick-film carbon electrodes** (ED-S1PE-C) are fabricated by **printing technologies** on a flexible and high-resist PET or a rigid ceramic substrate. These low-cost and disposable electrochemical sensors enable the use of **small sample volume**.

## » Thick-film based-electrode features

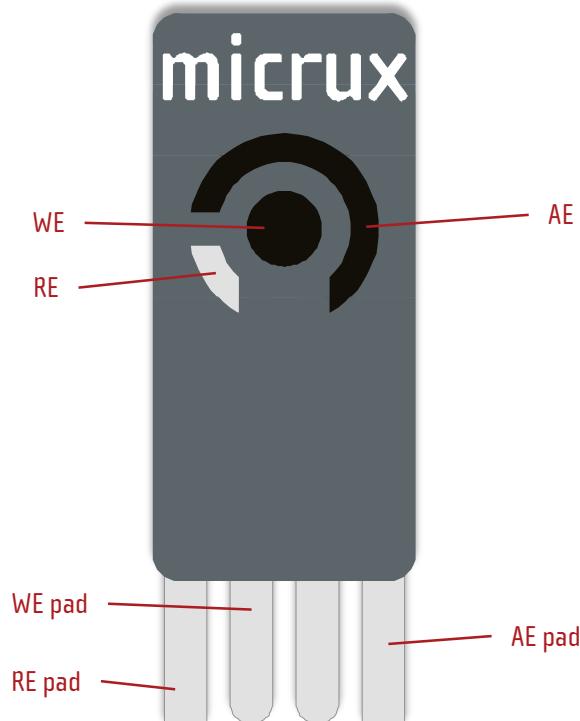

Printing technologies enable the manufacture of planar electrodes suitable for working with sample microdrop.

» <b>Standard dimensions:</b>	27.5 x 10.1 mm
» <b>Substrate:</b>	PET (white) / Ceramic ( $\text{Al}_2\text{O}_3$ )
» <b>Substrate thickness:</b>	350 $\mu\text{m}$ / 380 $\mu\text{m}$
» <b>WE dimensions:</b>	3 mm Ø (7,1 $\text{mm}^2$ )
» <b>Sample volume:</b>	20 – 50 $\mu\text{L}$
» <b>Electrode material</b>	
<b>Working electrode (WE):</b>	Carbon
<b>Reference electrode (RE):</b>	Silver (or Ag/AgCl)
<b>Auxiliary electrode (AE):</b>	Carbon

## » Thick-film electrode packs

Thick-film S1PE electrodes are supplied in **50 units packs**. They should be stored at room temperature in a dry place.

## » Applications


**Printed electrodes** are a suitable tool for **multiple applications**, providing many advantages such as low-cost, disposable, low reagent consumption as well as non-tedious pre-cleaning procedures.

Electroanalysis	Nanotechnology	Biosensors	Flow Analysis Systems
<ul style="list-style-type: none"> <li>✓ Study EC reactions</li> <li>✓ Trace EC analysis</li> <li>✓ In-vivo measurements</li> </ul>	<ul style="list-style-type: none"> <li>✓ Modified electrodes</li> <li>✓ New nanostructures</li> <li>✓ New nanomaterials</li> </ul>	<ul style="list-style-type: none"> <li>✓ EC transducers</li> <li>✓ New recognition elements</li> <li>✓ POC / wearable systems</li> </ul>	<ul style="list-style-type: none"> <li>✓ FIA Systems</li> <li>✓ HPLC</li> <li>✓ Capillary Electrophoresis</li> </ul>



## » Electrochemical cell

Carbon ([Ref. ED-51PE-C](#)) thick-film electrochemical sensors are based on a classical three-electrodes (working – WE, reference – RE and auxiliary – AE) approach.

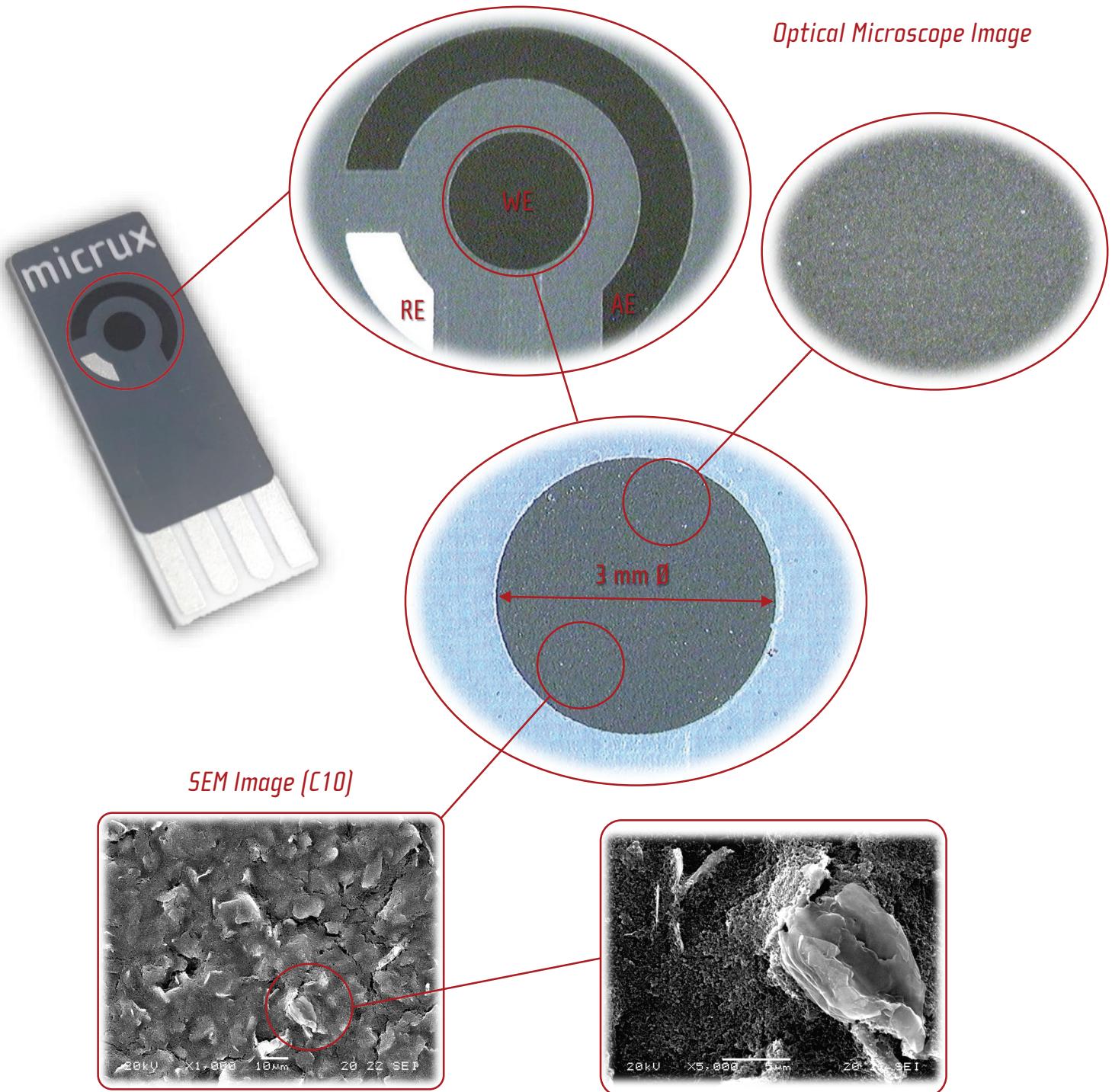


**Printed-based** electrodes are very useful in order to avoid tedious polishing of traditional solid electrodes, and make easier the development of **chemical-sensors** and **bio-sensors** for field analysis.




## » Thick-film carbon electrodes performance

» PRECISION INTRA-ELECTRODE

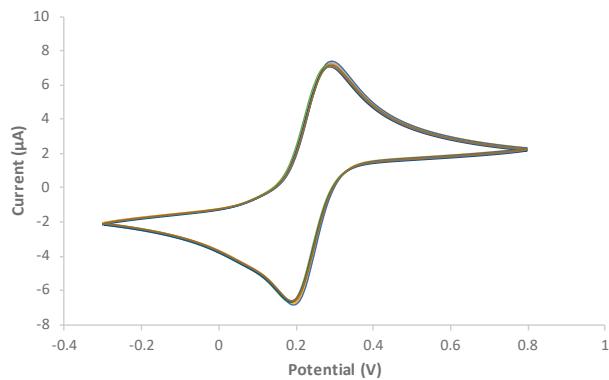


Successive cyclic voltammograms for 1 mM  $K_4Fe(CN)_6$  in 0.1 M  $H_2SO_4$  at the **same** thick-film carbon electrode [ED-51PE-C10].  $v = 50 \text{ mV/s}$ ,  $n = 10$ , RSD = 2%

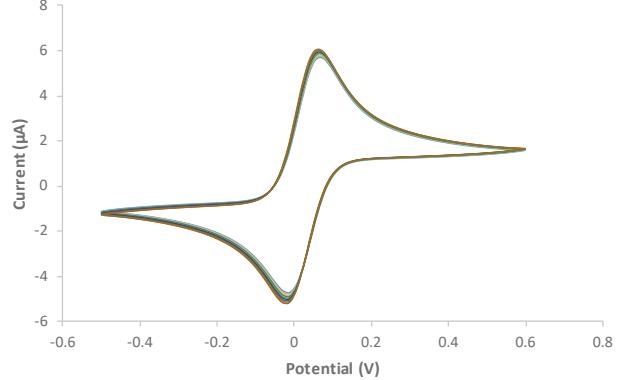


Successive cyclic voltammograms for 1 mM ferrocene methanol in 0.1 M  $H_2SO_4$  at the **same** thick-film carbon electrode [ED-51PE-C10].  $v = 50 \text{ mV/s}$ ,  $n = 10$ , RSD = 2%

» PRECISION INTER-ELECTRODE

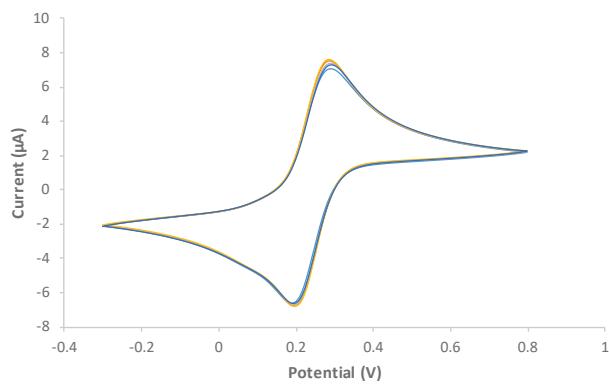


Cyclic voltammograms for 1 mM  $K_4Fe(CN)_6$  in 0.1 M  $H_2SO_4$  at **different** thick-film carbon electrodes [ED-51PE-C10].  $v = 50 \text{ mV/s}$ ,  $n = 5$ , RSD = 3%

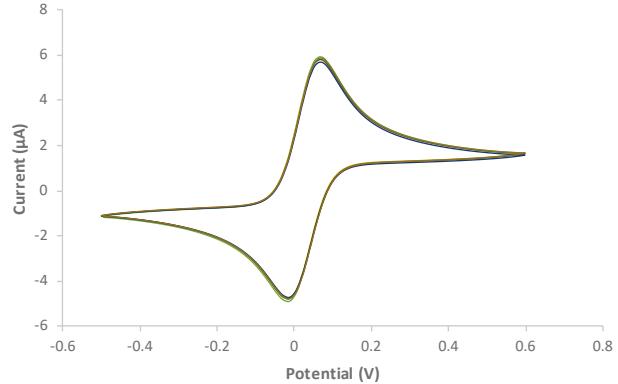


Cyclic voltammograms for 1 mM ferrocene methanol in 0.1 M  $H_2SO_4$  at **different** thick-film carbon electrodes [ED-51PE-C10].  $v = 50 \text{ mV/s}$ ,  $n = 5$ , RSD = 2%

» SWEEP RATE

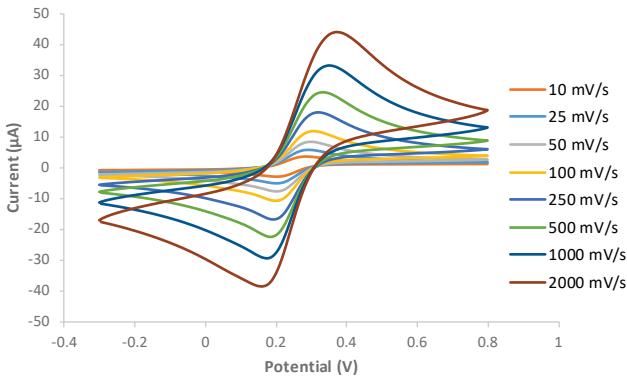

Cyclic voltammograms for 1 mM  $K_4Fe(CN)_6$  in 0.1 M  $H_2SO_4$  using different **scan rates** at a thick-film carbon electrode.

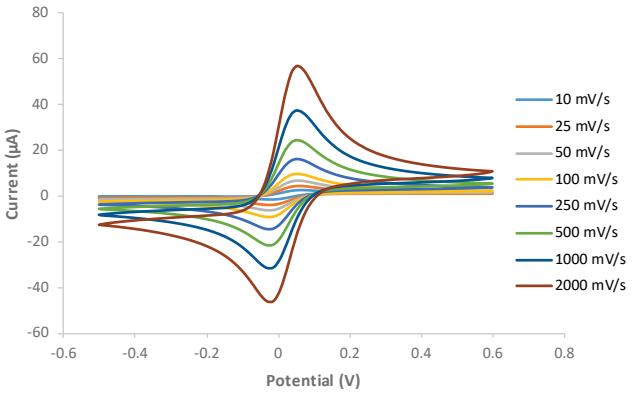


Cyclic voltammograms for 1 mM ferrocene methanol in 0.1 M  $H_2SO_4$  using different **scan rates** at a thick-film carbon electrode.

## » Thick-film electrodes selection

Screen-printed single electrodes (S1PE) are available in different materials (substrates and electrodes).

SKU	Substrate	WE	RE	AE
» ED-S1PE-C10	PET	Carbon	Silver	Carbon
» ED-S1PE-C10C	Ceramic	Carbon	Silver	Carbon
» ED-S1PE-C20	PET	Carbon	Silver	Carbon
» ED-S1PE-C20C	Ceramic	Carbon	Silver	Carbon
» ED-S1PE-C21	PET	Carbon	Silver/Silver Chloride	Carbon
» ED-S1PE-C21C	Ceramic	Carbon	Silver/Silver Chloride	Carbon
» ED-S1PE-CG20	PET	Graphene	Silver	Graphene
» ED-S1PE-CG21	PET	Graphene	Silver/Silver Chloride	Graphene
» ED-S1PE-Au10	PET	Gold	Silver	Gold
» ED-S1PE-Au10C	Ceramic	Gold	Silver	Gold

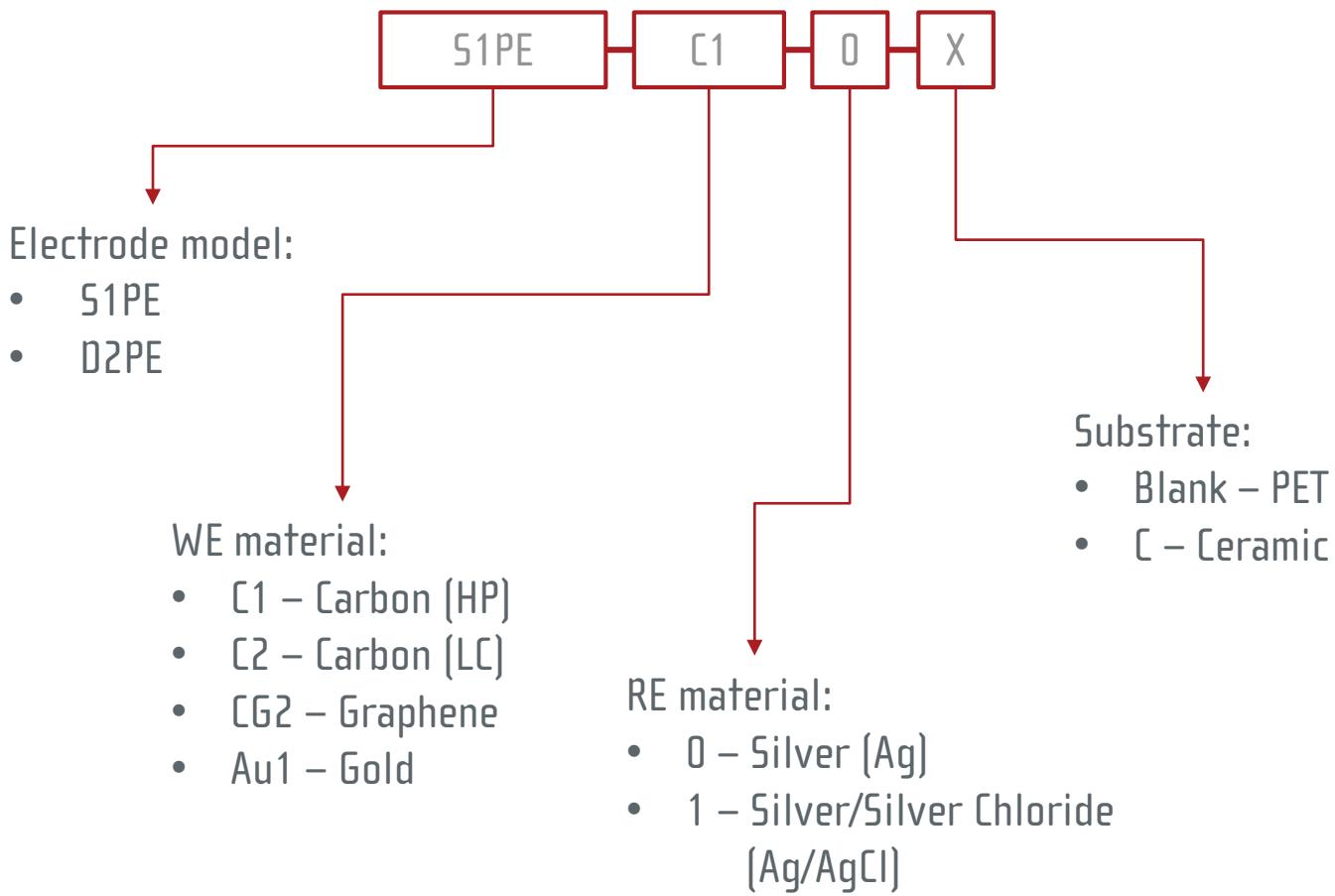

*Note: All combinations are NOT available. Please always check the availability for your specific needs.*



# Thick-film electrodes



## » Thick-film electrodes related accessories

Different **connectors** for interfacing the printed electrodes with any commercial potentiostat are also available at MicruX.

### » CABLE connector (ED-SPE-CABLE)

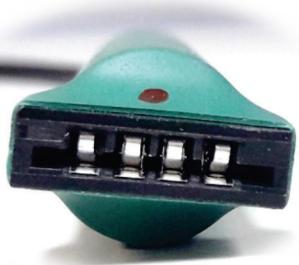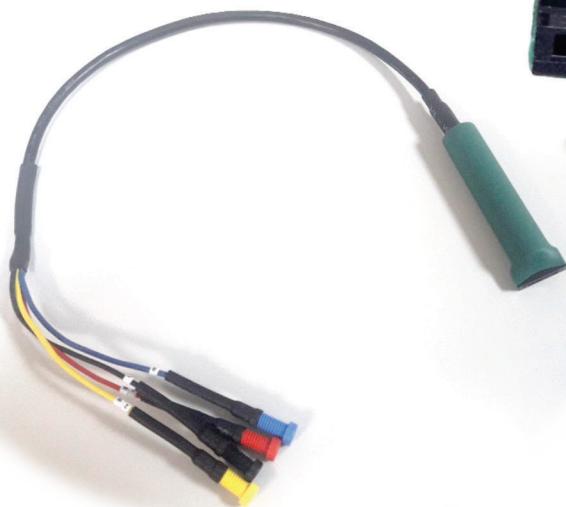

The **CABLE connector** (*Ref. ED-SPE-CABLE*) provides an interface between the electrodes (up to four contact pads) with the potentiostat, enabling the use of microvolume (20 – 50 µL sample drops) or dipping into a solution. The cable ends are available with **2 mm female or male bananas**.

*Dimensions: 50 cm long*

### » BOX Connector (ED-SPE-BOX)



The small **BOX connector** (*Ref. ED-SPE-BOX*) provides an interface between the electrodes (up to four contact pads) with any kind of potentiostat, enabling the use of microvolume (20 – 50 µL sample drops). The interface ends are available with **2 mm female bananas**.

*Dimensions: L58 x W40 x H15 mm*



Mora-Garay Industrial Park  
Juan de la Cierva, 2C, Bldg. # 6  
33211 · Gijón (Asturias) · SPAIN

Phone/FAX: +34 984151019

E-mail: [info@micruxfluidic.com](mailto:info@micruxfluidic.com)  
Web: [www.micruxfluidic.com](http://www.micruxfluidic.com)

